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GB Instruction Manual

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CAUTION!

Sharp-edged and pointed tools are often used when working with this device. Therefore, keep this device and all accessories and tools in a place that is inaccessible to children.

**Only allow children to work with the device under supervision!
Keep packaging material (plastic bags, rubber bands, etc.) away from children!**

Delivery content (Fig. 1)

Microscope (A), 2x/4x Objective (pre-assembled) (B), Power adapter with power cable (C), 2 Object plates (1x glass, 1x black/white) (D), 1 pair rubber eyecups (E), 1 pair eyepieces (WF 10x) (F), 1 pair eyepieces (WF 20x) (G), hook key (H), dust cover (I)

All parts (Fig. 2-6)

- 1 Eyepieces
- 2 Clamping screw (eyepiece)
- 3 Dioptre setting
- 4 Binocular head
- 5 Knurled screws (binocular head)
- 6 Focus wheel
- 7 Objektive holder
- 8 Objektive
- 9 Illumination (Incident/transmitted light)
- 10 Objecte clamps
- 11 On/Off switch
- 12 Dimmer wheel
- 13 Object plate
- 14 Selector swtch (illumination type)
- 15 Microscope head
- 16 Boreholes (Friktion setting)
- 17 Power socket
- 18 Battery compartment
- 19 Hook key (Friktion seetting)

1. Choosing a place for the microscope

Before starting, choose a sufficient location for microscoping. Please take care that your microscope stands on an even and solid ground. A power supply socket is required to charge the batteries.

Position your device so that it can be disconnected from the power supply at any time. The wall socket should always be located near the device and be easily accessible, since the plug on the power cord serves as a disconnecting device for the power supply.

2. Putting into operation**2.1 Power supply**

This microscope can be operated via mains power (230V/50Hz) or with batteries.

NOTE:

The microscope automatically detects the type of power supply. If the 230V mains power supply is available, this will be used. If only batteries with sufficient capacity are inserted, the microscope switches to battery operation.

2.1.1 Operation using mains power supply

1. Insert the barrel plug of the power adapter (C) into the power connection socket (18) on the back of the microscope.
2. Insert the Euro plug of the power adapter (C) into a suitable mains power socket (230V/50Hz).
3. Move the On/Off switch (11) to position ‚I‘ enable the power supply.
4. Move the On/Off switch (11) to position ‚0‘ disable the power supply.

2.1.2 Operation using batteries

1. Remove the cover of the battery compartment (13) on the bottom of the microscope.

2. Insert 3 AA/LR6-type batteries and replace the cover.

IMPORTANT:

When inserting the batteries, make sure that the battery poles are correctly aligned as indicated in the battery compartment!

3. Move the On/Off switch (11) to position ,I' enable the power supply.

4. Move the On/Off switch (11) to position ,0' disable the power supply.

NOTE:

This microscope does not have a charging function. Therefore, NEVER use any rechargeable batteries (accumulators)! Only use AA/LR6-type batteries.

2.2 Setting the illumination

This microscope can be operated with different types of LED illumination (incident light, transmitted light, combined).

1. Turn the selector switch (14) to the position for the desired illumination type: I = incident light, II = transmitted light, III = incident and transmitted light combined.

2. Turn the dimmer wheel (12) to adjust the brightness of the illumination.

3. Turn the selector switch (14) to the '0' position to switch the illumination off completely.

NOTE:

The LED illumination has a service life of approximately 20,000 operating hours. If a defect occurs, please contact the service center.

3. Observation

First, remove the dust protective cover (6) from the objective holder (5). Place the desired object in the middle of the stage (8) and secure it with the clamps (9) if necessary. The objective's magnification (2x or 4x) can be adjusted by turning the objective holder (5). It is recommended to start observations always with the lowest magnification. Now look through the eyepieces (1) at the binocular head (3). Adjust the sharpness with the focuser (4).

4. Focusing and dioptre compensation

Set the dioptre scale (2) to „0“ and observe an object. Now close your left eye and adjust the focus with the focuser (4) for your right eye. Then, close your right eye and adjust the focus for your left eye with the dioptre scale (2). Now, the microscope is adjusted for your eyes. Note the setting of the dioptre scale you need to use in future.

5. Eye distance

The Researcher ICD is equipped with an adjustable binocular head. To set your eye distance, move the eyepieces against each other until you found the ideal position.

6. Replacing the Eyepieces

To replace the eyepieces you need to loosen the clamp screw on the underside of the eyepiece holder (17). You can then easily remove and replace the eyepieces.

The clamp screws serve only as protection so that the eyepieces will not fall out.

NOTE:

The supplied rubber eyecups (E) can be placed over the eyepieces to protect against stray light from the side. This also provides a more comfortable viewing experience for the user.

7. Adjusting the friction of the focusing drive

It can happen that the Microscope head is no longer held by the focusing drive and shuts down because of his net weight.

With the BRESSER Researcher ICD you can adjust the friction with the enclosed wrench (19).

The wrench is simply hooked into two holes of the wider shaft (18). By rotating you can adjust the friction strongly or lightly.

8. Cleaning and maintenance

Your microscope is a high-quality optical instrument. Please take care that neither dust nor fingerprints take place on the optical surfaces. If necessary, clean the microscope body with a slightly moistened soft cloth. Use a separate clean cloth with some drops of alcohol to clean the optical surfaces. After using, the microscope should be protected with its dust cover.

NOTE:

Please consider that a sufficient treated microscope keeps its value and optical quality for many years.

9. Technical data

Binocular head

Eyepieces: 10x and 20x (2 pcs. each), 30.5 mm diameter

Objectives: 2x and 4x, Stereo

Total magnification: 20x, 40x, 80x (The total magnification can be calculated by multiplying the objective magnification with the eyepiece magnification, e.g.. $2 \times 40 = 80$)

Lighting: LED

Power supply: 230V/50Hz mains power via supplied power adapter or 3/6 AA/LR6 type batteries

11. CE Declaration of Conformity

 Bresser GmbH has formed a „Conformity Declaration“ in line with the applicable directive. This declaration is available for review under www.bresser.de/download/5803100/CE/5803100_CE.pdf

12. Disposal

Dispose of the packaging materials separately.

When disposing of the device, observe the current legal regulations!

Information on proper disposal is available from the municipal waste disposal service providers or the environmental agency..

Do not dispose of electric equipment together with household waste material!

In observance of European Directive 2002/96/EC on waste electrical and electronic equipment (WEEE) and its implementation in accordance with national law, electric equipment that have reached the end of its life must be collected separately and returned to an environmentally compatible recycling facility.

13. Warranty & Service

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full warranty terms as well as information on extending the warranty period and details of our services at www.bresser.de/warranty_terms.

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GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

E-Mail: service@bresseruk.com
Telephone*: +44 1342 837 098

BRESSER UK Ltd.
Suite 3G, Eden House
Enterprise Way
Edenbridge, Kent TN8 6HF
United Kingdom

*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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Bresser UK Ltd.