## **Biolux NV**

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*Expand your horizon*

## General Information

We wish you much success in your work with your new BRESSER Biolux NV microscope.

### About this Instruction Manual

Please read the safety instructions in this manual carefully. To avoid damage to the unit and the risk of injury, only use this product as described in the manual.

Keep the instruction manual handy so that you can easily look up information on all the functions.

#### **DANGER!**

You will find this symbol in front of every section of text which deals with the risk of severe injury or even death in the event of improper use.

You will find this symbol before every section of text that deals with the risk of minor to severe injuries resulting from improper use.

#### **NOTE!**

You will find this symbol in front of every section of text which deals with the risk of damage to property or the environment.

### Intended Use

This product is intended only for private use.

It was developed for the magnified display of things in nature.

## General Warning

#### **DANGER!**

Tools with sharp edges and points are often used when working with this device. For this reason, store this device and all accessories and tools in a location that is out of the reach of children. There is a RISK OF INJURY!

#### **DANGER!**

This device contains electronic components which operate via a power source (batteries). Do not leave children unattended when using the device. Only use the device as described in the manual, otherwise you run the RISK of an ELECTRIC SHOCK.

#### **DANGER!**

Do not expose the device to high temperatures. Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.

Keep batteries out of the reach of children! Make sure you insert the batteries correctly. Empty or damaged batteries could cause burns if they come into contact with the skin. If necessary, wear adequate gloves for protection.

#### **DANGER!**

Use only the recommended batteries. Always replace weak or empty batteries with a new, complete set of batteries at full capacity. Do not use batteries from different brands or with different capacities. The batteries should be removed from the unit if it has not been used for a long time.

The manufacturer is not liable for damage related to improperly installed batteries.

#### **DANGER!**

Children should only use the device under adult supervision. Keep packaging material (plastic bags, rubber bands, etc.) out of the reach of children! There is a RISK OF CHOKING!

**CAREFUL!**

Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after use. In the event of contact with the eyes or mouth, rinse thoroughly with water. In the event of pain, contact a doctor immediately and take the substances with you.

**NOTE!**

Do not disassemble the device. In the event of a defect, please contact your dealer. He will contact the Service Centre and can send the device in to be repaired, if necessary.

## Operating instructions

### Parts overview (Fig. 1+2):

- ❶ 5x WF Eyepiece
- ❷ 16x WF Eyepiece
- ❸ Barlow lens
- ❹ MicrOcular
- ❺ Eyepiece holder
- ❻ Microscope head
- ❼ Objective-revolver
- ❽ Microscope stage
- ❾ Focus wheel
- ❿ LED lighting (transmitted light)
- ⓫ Microscope housing
- ⓬ Microscope base
- ⓭ Battery compartment (3x AA)
- ⓮ 5 Slides, 10 Covering glasses and 5 preparations in a plastic box
- ⓯ Dimmer
- ⓰ Colour filter disc
- ⓱ LED lighting (direct light)
- ⓲ Direct light / transmitted light switch
- ⓳ Microscope tools:
  - a) Pipette; b) Tweezers
- ⓴ Prawn breeding plant
- ⓵ MicroCut
- ⓶ Specimens:
  - a) yeast; b) Gum media (specimen inclusion medium); c) sea salt;
  - d) Prawn eggs
- ⓷ Carrying case
- ⓸ Locking screw
- ⓹ cross stage

### 1. General/Location

First you must make sure that your microscope is on a stable and solid surface.

### 2. Electric LED lighting with dimmer

Check before use whether the light switch (Fig. 1,18) is set to "off".

The microscope has two lighting units. Lighting can be of three types. Set the switch (Fig 1, 18) to „II“ to light the specimen from above (reflected light) or „I“ to light it from below (direct light). Use setting „III“ to have the specimen simultaneously subjected to direct and transmitted light. The transmitted light unit (Fig 1, 10) is used for transparent specimens (those on glass slides). To view solid non-transparent specimens use the direct light unit (Fig 1, 17). Use of both forms of lighting simultaneously is only recommended for semi-transparent specimens. This operating mode is not recommended for direct light specimens on slides as it may cause reflection on the slide.

Use the switch (Fig 1, 18) to select the desired lighting mode and set the dimmer to the desired brightness (Fig. 1, 15).

As your device has continuously controllable lighting (dimmer) optimal illumination of the object to be viewed is guaranteed.

### 3. Colour filter disc

The colour filter under the microscope stage (Fig. 1, 8) aids in viewing very bright and transparent objects. Just select the right colour for the specimen in question. The components of colourless/transparent objects (e.g. starch particles, single-cell -specimens) can thus be better recognised.

### 4. Attach cross stage

A stage is included along with your microscope (located in the case). To install this on the microscope, proceed as follows:

1. Place the cross stage (Fig. 4, 25) on the microscope stage (Fig. 4, 8) in such a way that the attachment screws (Fig. 4, D) are positioned over the middle threaded hole (Fig. 4, E).
2. Tighten the attachment screws securely.

### 5. Microscope setup

The microscope's eyepiece (Fig 1, 6) will now be prepared for the first observation.

First, loosen the locking screw (Fig. 1, 24) and rotate the eyepiece into a convenient position.

Begin every observation with the lowest magnification. Place the microscope's stage (Fig. 1, 8) with the focus knob (Fig 1, 9) into the lowest position and rotate the objective revolver (Fig. 1, 7) until it locks on the lowest magnification (4x).

#### NOTE:

First, drive the Microscope's stage (Fig. 1, 8) in its lowest position before changing the objective in order to prevent damages.

Insert the 5x eyepiece (Fig. 3, 1) in the Barlow lens (Fig. 3, 3).

Take care, that the Barlow lens is inserted completely in the monocular head (Fig. 3, 5).

### 6. Observation

After you have set up the microscope with the corresponding illumination, the following principles are important:

Begin each observation with a simple observation at lowest magnification, so that the centre and position of the object to be viewed is in focus.

The higher the magnification the more light is required for good picture quality.

Place a permanent slide culture (Fig. 5, 14) directly under the microscope lens on the plate (Fig. 5, 8) and clamp it on the cross stage (Fig. 5, 25). To do so push the lever (Fig. 5, C) aside. The specimen to be examined must be directly over the lighting. If not turn the two knurled screws (Fig. 5, A+B) on the cross stage.

#### TIP:

On the cross stage (Fig 5, 25) there are two knurled screws (Figs 5, A+B). They are used to precisely position the specimen laterally (Fig 5, A) and vertically (Fig 5, B).

Look through the eyepiece (Fig. 1, 1/2) and turn carefully the focusing wheel (Fig. 1, 9) until you can see a sharp picture.

Now you can get a higher magnification, while you pull out slowly the Barlow lens (Fig. 6, 3) of the monocular barrel (Fig. 6, 5). With nearly entirely pulled out Barlow lens the magnification is raised to 2x.

For still higher magnification you can put the 16x eyepiece (Fig. 2, 2) into the objective revolver (Fig. 1, 7) and set on higher position (10x / 40x).

#### TIP:

Depending on the preparation higher magnifications do not always lead to better pictures.

With changing magnification (eyepiece or objective lens changes, pulling out of the Barlow lens) the sharpness of the image must be newly defined by turning the focusing wheel (Fig. 1, 9).

**NOTE:**

Please be very careful when doing this. When you move the microscope stage upwards to fast the objective lens and the slide can touch and become damaged.

## 7. Viewed Object – condition and preparation

### 7.1. Condition

With the Barlow lens nearly fully extended magnification can be doubled. Both transparent and non-transparent specimens can be examined with this microscope, which is a direct as well as transmitted light model. If opaque specimens are examined - such as small animals, plant parts, tissue, stone and so on - the light is reflected from the specimen through the lens and eyepiece, where it is magnified, to the eye (reflected light principle, switch position I). If opaque specimens are examined the light from below goes through the specimen, lens and eyepiece to the eye and is magnified en route (direct light principle, switch position II). Many small organisms of the water, plant parts and finest animal components have now from nature these transparent characteristic, other ones must be accordingly prepared. Is it that we make it by means of a pre-treatment or penetration with suitable materials (media) transparent or thus

that we cut finest wafers off of them (hand cut, MicroCut) and these then examine. With these methods will us the following part make familiar.

### 7.2. Creation of thin preparation cuts

Specimens should be sliced as thin as possible, as stated before. A little wax or paraffin is needed to achieve the best results. A candle can be used for the purpose. The wax is put in a bowl and heated over a flame. The specimen is then dipped several times in the liquid wax. The wax is finally allowed to harden. Use a MicroCut (Fig 2, 21) or knife/scalpel (carefully) to make very thin slices of the object in its wax casing.

**DANGER!**

Be extremely careful when using the knives/scalpels or the MicroCut. There is an increased risk of injury due to the sharp edges!

These slices are then laid on a glass slide and covered with another.

### 7.3. Creation of an own preparation

Put the object which shall be observed on a glass slide and give with a pipette (Fig. 8, 19a) a drop of distilled water on the object (Fig. 8).

Set a cover glass (in each well sorted hobby shop available) perpendicularly at the edge of the water drop, so that the water runs along the cover glass edge (Fig. 9). Lower now the cover glass slowly over the water drop.

**TIP:**

The gum medium supplied (Fig 2, 22b) is used to make permanent slide cultures. Add it instead of distilled water. The gum medium hardens so that the specimen is permanently affixed to its slide.

## 8. Experiments

If you made yourself familiar with the microscope already, you can accomplish the following experiments and observe the results under your microscope.

### 8.1. Newspaper print

#### *Objects:*

1. A small piece of paper from a newspaper with parts of a picture and some letters
2. A similar piece of paper from an illustrated magazine

Use your microscope at the lowest magnification and use the preparation of the daily paper. The letters seen are broken out, because the newspaper is printed on raw, inferior paper. Letters of the magazines appear smoother and more complete. The picture of the daily paper consists of many small points, which appear somewhat dirty. The pixels (raster points) of the magazine appear sharply.

### 8.2. Textile fibres

#### *Objects and accessories:*

1. Threads of different textiles: Cotton, line, wool, silk, Celanese, nylon etc..
2. Two needles

Each thread is put on a glass slide and frayed with the help of the two needles. The threads are dampened and covered with a cover glass. The microscope is adjusted to a low magnifica-

tion. Cotton staples are of vegetable origin and look under the microscope like a flat, turned volume. The fibres are thicker and rounder at the edges than in the centre. Cotton staples consist primary of long, collapsed tubes. Linen fibres are also vegetable origin; they are round and run in straight lines direction. The fibres shine like silk and exhibit countless swelling at the fibre pipe. Silk is animal origin and consists of solid fibres of smaller diameter contrary to the hollow vegetable fibres. Each fibre is smooth and even moderate and has the appearance of a small glass rod. Wool fibres are also animal origin; the surface consists of overlapping cases, which appear broken and wavy. If it is possible, compare wool fibres of different weaving mills. Consider thereby the different appearance of the fibres. Experts can determine from it the country of origin of wool. Celanese is like already the name says, artificially manufactured by a long chemical process. All fibres show hard, dark lines on the smooth, shining surface. The fibres ripple themselves/crinkle after drying in the same condition. Observe the thing in common and differences.

### 8.3. Salt water prawns

#### 8.3.1 Accessories

1. Shallow vessel for approx. 2-3 L (canning jar, glass bowl, mini aquarium or Faunabox) - to ensure sufficient air exchange at the water surface.

2. 2-3 liters of rainwater, Volvic (or other still water)
  3. Prawn eggs (Fig. 2, 22d)
  4. Sea salt (preferably uncleaned, without additives, about 100g)
- Note:** the supplied amount of sea salt (Fig. 2, 22c) is not sufficient for this long-term approach!
5. Yeast (Fig. 2, 22a) – Food
  6. Pipette (Fig. 2, 19a) or a thin tube - to remove animals or change water
  7. Prawn breeding plant (Fig. 2, 20) – to observe the animals under the microscope or to let only a few nauplii hatch (not recommended as it is not possible to keep them in this vessel)
  8. Optional: sodium bicarbonate "natron" (approx. 5g)

#### 8.3.2. The lifecycle of the saltwater prawn

The saltwater prawn or artimia salina has an unusual and interesting life cycle for scientists. The female's eggs are hatched without the need for male crabs to fertilize them. The resulting baby crabs are all female. Under unusual circumstances, such as when a marsh dries up, the eggs may produce male crayfish. These males fertilize the eggs of the females, resulting in a specific type of egg. These eggs, called "permanent eggs" or cysts, have a thick shell that protects the egg. The "winter eggs" are very resilient and remain viable even if the marsh or lake dries up, causing the death of the entire crayfish population. They can

remain in a "dormant" state for 5-10 years. The eggs will hatch as soon as the proper environmental conditions return. The supplied crayfish eggs are of this nature.

### 8.3.3. Hatching of the salt water prawn

To hatch, raise and observe the crayfish for a long time, it is important to first prepare a large batch of saline solution that corresponds to the living conditions of the crayfish. For this purpose you need a vessel for about 2-3 L (canning jar, glass bowl, mini aquarium or Faunabox). For the preparation it is best to use a rather shallow vessel to ensure sufficient air exchange at the water surface.

The permanent eggs are prepared with 2-3 liters of rainwater. Add one tablespoon (about 10-15 g) of sea salt without additives per 0.5 liters of water. Unpurified, gray sea salt is optimal. To make the pH value slightly alkaline (pH about 7.5-8.0), a pinch of sodium hydrogen carbonate "natron" can be added per 0.5 L. The sodium hydrogen carbonate is then added to the water. If the water is chlorinated or contains copper (e.g. from pipes, flow heater), you can use mineral water low in minerals and pH neutral (e.g. Volvic) for the preparation, since invertebrates react very sensitively to chlorine and heavy metals. The vessel should be kept at room temperature in a bright place, but not directly in the sun. After 24-48 hours at room temperature,

the first 0.3 mm to 0.5 mm larvae, called nauplii, hatch. These paddle through the water with the rudder antennae. Initially, the Artemia moult at very short intervals (a few hours). Only after several molts does the number of leg pairs increase. Depending on the ambient temperature, the larvae are fully grown in 6 to 10 weeks. They will soon have bred a whole generation of saltwater crayfish that are constantly reproducing.

### 8.3.4. The saltwater prawn under the microscope

The animals can be optimally observed with a reflected light microscope. To do this, place a few animals in some water in the saltwater crayfish incubator. Smaller animals can also be pipetted with a pipette with the largest possible opening or aspirated with a thin tube.

You can remove the lid of the incubator and view it in its entirety under the microscope.

During microscopy, the specimen may heat up due to the microscope illumination. Important: Keep observation times short. Switch off the illumination during breaks and remove the brine shrimp incubator with the animals from the microscope! After observation, quickly return the animals to the culture vessel.

### 8.3.5. Feeding and water change

To keep the saltwater prawn alive, they must be fed occasionally. This should be done carefully,

otherwise the water will spoil and the crayfish will be poisoned. Feeding is best done with dry yeast powder (Fig. 2, 22a), spirulina powder or "green water". Give some to the crayfish starting on the 2nd day after hatching and then every other day. You can, in order to preserve the culture as long as possible, siphon off accumulating mulm and food residues on the bottom with a pipette. Fill up the evaporated water with deionized or distilled water. Every two weeks you should also replace a part of the water. It is important that the water change is done carefully; the fresh water should have the same temperature and composition, and be added over a period of at least 1-2 hours. With a little luck, this way the culture can be maintained for a period of up to 6 months (max. age). On the soil, with a little luck, you can find new cysts that can be dried. From them you can hatch a new generation.

**ATTENTION!**  
SALTWATER PRAWN AND  
THEIR EGGS ARE NOT  
SUITABLE FOR CONSUMPTION!

For more information on primeval crabs, visit our website at:

<https://www.bresser.de/c/en/support/ratgeber/mikroskopie/urzeitkrebse/>

## 9. Use of the camera with suitable Software

### 9.1. Install Bresser CamLab Lite software once

#### *Free download for the latest Bresser CamLab Lite software*

The free download for the current Bresser CamLab Lite software version can be found on our product page at:

<https://www.bresser.de/P5116200>

Then select the "Software" tab. Here you will find the download links for the operating systems Windows and MacOS. You will also find the software download information in printed form in the case (Fig. 2/23).

#### *CamLab Lite software tutorial*

The operating instructions for the CamLab Lite software can be found on our product page at: [www.bresser.de/P5116200](http://www.bresser.de/P5116200)

In the "Instructions" tab, the operating instructions are available as a pdf. In the "Instructions" tab, the operating instructions are available as a pdf.

### 9.2. Using the ocular camera for observations on the PC

#### **ATTENTION!**

BEFORE using the micrOcular, make all settings on the microscope according to section 6. Make sure to place a specimen with as much contrast as possible on the microscope (e.g. stained plant cuttings).

#### **NOTE:**

The MicrOcular only works without the supplied Barlow lens! The magnification setting is changed by using the MicrOcular and must be readjusted using the focus wheel.

Remove the Barlow lens (Fig. 6, 3) and the eyepiece currently in use from the eyepiece tube (Fig. 6, 5). Instead, insert the micro eyepiece (Fig. 7, 4) into the eyepiece tube (Fig. 7, 5) as shown in Fig. 7.

### 9.3. Photo and video recording via CamLab Lite software

Connect the micrOcular to your computer. To do this, use the appropriate cable to connect the camera module to a USB port.

Then open the previously installed CamLab Lite software. The software initially opens without displaying the live view of the camera. At the top of the left side you will find the first menu item "Camera list". Click with the left mouse button on the name of the camera MikrOkularHD. Now the live view of your high-contrast specimen opens. Refocus the specimen to obtain a sharp image in the preview. The entire image field is always displayed in the software.

#### **NOTE:**

The camera shows a smaller section than can be seen in the eyepiece. This is normal. If necessary, select a different magnification on the microscope to capture a suitable image section.

The recording is started via the "Snapshot" (single image) or "Record" (video) button. The storage location and the language of the software can be selected via the setting wheel. Select the resolution of the preview or the recording via the "Live" menu. The image brightness is automatically controlled. Make sure that the microscope illumination is set as bright as possible. If you would like to control the image brightness man-

ually, remove the check mark from "Auto Exposure" and operate the slider at "Exposure Time".

**NOTE:**  
For an optimal result, settings can also be made in the "Color Balance" menu. Depending on the camera model, individual controls may be inactive if the camera does not support these functions.

**TIP:**  
To work permanently with the micro eyepiece, it is recommended to always operate the micro eyepiece on one and the same USB port.

### 10. Care and maintenance

Disconnect the device from the power source before cleaning (remove batteries)! Clean the device only externally with a dry cloth.

**NOTE:**  
Do not use cleaning fluid to avoid damage to the electronics.

Only clean the lenses (eyepieces and/or objectives) with a soft and lint-free cloth (e.g. microfiber).

**NOTE:**  
Do not press the cloth too hard to avoid scratching the lenses.

To remove heavier dirt residues, moisten the cleaning cloth with an eyeglass cleaning liquid and wipe the lenses with it using little pressure.

Protect the device from dust and moisture! Allow it to acclimatize at room temperature for some time after use - especially in the case of high humidity - so that the residual moisture can be reduced.

### 11. Troubleshooting

Error	Solution
<b>No image recognizable (when observed by eye)</b>	<ul style="list-style-type: none"> <li>• Switch on light</li> <li>• Check that the selected objective is engaged in its end position on the revolving nosepiece</li> <li>• Readjust the image sharpness</li> <li>• Check if the specimen is placed in the center of the image field</li> <li>• Check the brightness setting</li> <li>• Check whether the illumination setting on the incident/transmitted light selector switch has been selected appropriately</li> </ul>
<b>Image is shaded sideways</b>	<ul style="list-style-type: none"> <li>• Check whether the color filter disk is locked in the desired end position</li> </ul>
<b>No live image recognizable in the camera software on the PC</b>	<ul style="list-style-type: none"> <li>• Under menu item "Camera list" click with the left mouse button on the name of the camera MikrOkularHD</li> </ul>

## 12. Specifications

### 12.1. System requirements

- **Windows 8.1, 10 or 11**
- **Intel i3, i5, i7 or i9 processor from 2nd generation onwards or AMD FX or Ryzen processor**
- **Optical Disk Drive or Internet connection for software download**
- **USB 2.0**

### 12.2 Magnification table

<i>Eyepiece</i>	<i>Objective</i>	<i>Magnification</i>	<i>with Barlow lens</i>
5x	4x	20x	40x
5x	10x	50x	100x
5x	40x	200x	400x
16x	4x	64x	320x
16x	10x	160x	320x
16x	40x	640x	1280x

## Declaration of Conformity

 Bresser GmbH has issued a „Declaration of Conformity“ in accordance with applicable guidelines and corresponding standards. This can be viewed any time under [https://www.bresser.de/download/5116200/CE/5116200\\_CE.pdf](https://www.bresser.de/download/5116200/CE/5116200_CE.pdf)

## UKCA Declaration of Conformity

 Bresser GmbH has issued a „Declaration of Conformity“ in accordance with applicable guidelines and corresponding standards. This can be viewed any time under [https://www.bresser.de/download/5116200/UKCA/5116200\\_UKCA.pdf](https://www.bresser.de/download/5116200/UKCA/5116200_UKCA.pdf)

## Disposal

 Dispose of the packaging materials properly, according to their type (paper, cardboard, etc). Contact your local waste disposal service or environmental authority for information on the proper disposal.

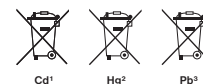
 Do not dispose of electronic devices in the household garbage!

As per the Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected sepa-

rately and recycled in an environmentally friendly manner.

 In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please make sure to dispose of your used batteries as required by law — at a local collection point or in the retail market. Disposal in domestic waste violates the Battery Directive.

Batteries that contain toxins are marked with a sign and a chemical symbol.

- <sup>1</sup> battery contains cadmium
- <sup>2</sup> battery contains mercury
- <sup>3</sup> battery contains lead

## Warranty & Service

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary warranty period as stated on the gift box, registration on our website is required.

You can consult the full warranty terms as well as information on extending the warranty period and details of our services at [www.bresser.de/warranty\\_terms](http://www.bresser.de/warranty_terms).

## Scope of delivery

- Solid metal Microscope
- mechanical desk
- HD camera with metal housing, resolution 1280×720
- 2 eyepieces: 5x and 16x
- 3 objectives: 4x, 10x and 40x
- 1 Barlow lens with 2x magnification
- Microscope tools
- cross stage
- 5 prepared slides
- 5 blank slides
- 3x AA batteries
- Dust cover and hard case

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### GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

E-Mail: [service@bresseruk.com](mailto:service@bresseruk.com)  
Telephon\*: +44 1342 837 098

#### **BRESSER UK Ltd.**

Suite 3G, Eden House  
Enterprise Way  
Edenbridge, Kent TN8 6HF  
Great Britain

\*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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