

# HDMI FULL HD DISPLAY (13.3") FOR MIKROCAM PRO HDMI AND MINI HDMI MANUAL Art. No. 5914120

**A** The connecting plate for connection to the microscope camera\* is located on the folding support of the display

Accessories:  
2x Mounting screws  
1x Allen (hey) key  
1x HDMI cable  
1x Y-cable hollow plug/USB-C

**B** The connection plate to the display is located on the side of the microscope camera\*

**C** Attach the display to the microscope camera\* and the microscope head\*

Fasten to the outside of the folding support using the two mounting screws

**D** Connect HDMI display via HDMI cable and establish power supply (use camera power supply\* (connection 3))

Left-sided interfaces

- 1a Not used
- 2x USB-C
- 2a HDMI

Right-sided interfaces (e.g. MikroCam MINI FULL HD HDMI):

- 2b HDMI
- 1b DC 12V

Connections:

- 2a: HDMI cable
- 2b: Y-cable hollow plug/USB-C
- 3: Y-cable hollow plug/USB-C

**E** Use the control buttons to adjust the display optimally to your needs

Right-sided interfaces and control buttons

- 4 Menu navigation
- 5 Long: Display ON/OFF  
Short: Menu (brightness, contrast)  
Not required (USB)

Manual/Accessories/Notes:  
[www.bresser.de/P5914120](http://www.bresser.de/P5914120)

\*Not included in the scope of delivery

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail:  
[service@bresseruk.com](mailto:service@bresseruk.com) · Telephone: +44 1342 837098  
BRESSER UK Ltd. · 3G Eden House · Enterprise Way · Edenbridge · Kent TN8 6HF · Great Britain